e-Loop

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Annealing without Flame!

The electrical sterilization system e-Loop is most suitable for the sterilization of inoculation loops, needles and instruments. It is ideal for laboratories and safety cabinets where the use of gases and open flames is not permitted.

Robust and temperature-shock resistant

The sterilizing tube, which is made of wear-resistant quartz glass, reaches its optimal temperature of 900°C (1650°F) after only a few minutes. To sterilize the inoculation loop, simply insert it into the quartz tube. After 5-7 seconds, the inoculation loop is sterilized. A spray protector is integrated for more safety when handling with pathogenic material. Even clogging is not an issue. Any residue remaining on the smooth surface of the quartz tube can be removed easily. A 100% stainless steel design makes the e-Loop resistant to even the most extreme laboratory conditions.*

Flexible and safe

The sterilization tube can lock itself into place, using the specially designed angular adjusting device. The low and stable housing facilitates ergonomic operation; the unique design protects the working surface from contamination. Exceptional passive safety feature: After prolonged use, the residual heat display signals a hot surface of the sterilization tube in order to protect the user from burns.

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e-Loop

with 7 angular adjustments
residual heat display
tray for 3 instruments

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- Stainless steel design
- Long - lasting sterilization tube (quartz glass)
- Sterilization period 5-7 seconds*
- Self - locking angular adjustment

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The range:

e-Loop
 with 7 angular adjustments
 Residual heat display
 Tray for 3 instruments, screwdriver
 for 230 V mains voltage **Art.-No. LF5.001.000**
 for 110 V mains voltage **Art.-No. LF5.002.000**

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Technical Data:

Diameter of quartz glass tube: Ø 15 mm
 Length of quartz glass tube: 118 mm
 Angular adjustments: 7
 Sterilization temperature: 900 °C - 950 °C (1650 °F - 1742 °F)
 Power supply: 120 W
 Heat transfer: a stainless steel surface, located above, warms up at motionless air and at an ambient temperature of 20 °C
 Distance: 20 cm: 55 °C / 30 cm: 38 °C / 40 cm: 32 °C
 Measurements (w x h x d): 175 x 129 x 139 mm
 Weight: 1100 g
 Licenses: CE: EN 61326-1, EN61000-3-2, EN 61010
 EEC guidelines: 89/336/EEC, 73/23/EEC

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Quartz glass steri. tube
 wear- persistent and
 temperature-shock
 resistant

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Support for all inoculation loop holders
 made of stainless steel (w/o Fig.)

Art.-Nr./Art.-No. LF5.000.100

Tray / Tray
 links montierbar,
 aus Edelstahl
 for left side,
 made of
 stainless steel

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Holding device exclusive for
Biotool inoculation loop holders
 made of stainless steel, suitable for all angles

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Inoculation loop holder
 made of stainless steel,
 with sleeve nut,
 for inoculation loops
 with wire Ø 0.6 - 1 mm

Art.-Nr./Art.-No.
LF5.000.360

73/23/EEC
 89/336/EEC

EN 61326-1 A1/A2/A3
 EN 61000-3-2:2000
 EN 61010

BIOTOOLSWISS

Biotool AG
 Mühlegasse 7
 CH - 3400 Burgdorf

Telefon: ++41 (0)34 423 50 60
 Email: info@biotoolswiss.com
 Internet: http://www.biotoolswiss.com