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Autoloop^{PRO}

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The range:

Autoloop^{PRO}

includes 4 inoculation loop holders
4 inoculation loops
Pedestal
Instruction manual
2-year warranty
Laboratory gas burner **not** included
Foot pedal connector on request

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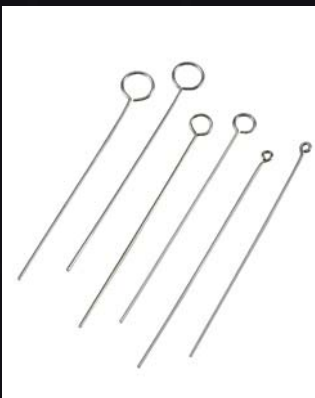
Spare inoculation loop holder
made of stainless steel,
with sleeve nut, for inoculation
loops with wire Ø 0.6 - 1 mm
Art.-Nr./Art.-No. 6.000.360

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Spare inoculation loops
made of special stainless steel,
wire Ø 0.6 mm

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Technical data:

Technology: Microprocessor
Display: fully graphic dot-matrix
Rotational direction: left / right, depends on removed holder
Inoculation loop holder: 1 - 4, with optical holder detection
Time of flame sterilizing: 1 - 15 sec
Time for cool-down period: 1 - 45 sec
Safety features: temperature control
flame activation warning
indicator of remaining cool-down
tilt sensor
residual heat display

Adjustable sterilization monitor: 1 - Off
2 - low intensity
3 - middle intensity
4 - high intensity
Stand by: automatic unit switch off: 1 h
(or with stand-by time of the burner)
Operating voltage: via Fuego (foot pedal socket)
Casing: stainless steel & aluminum,
UV and solvent resistant
Measurements: Pedestal (Art.-No. 8.000.400):
135 x 180 mm (w x d)
Pedestal (Art.-No. 6.000.400):
135 x 135 mm (w x d)
Height Autoloop^{PRO}: 305 mm
Weight: Art.-No. 8.000.400: 1200 g
Art.-No. 6.000.400: 1100 g
CE: EN 61326-1, EN61000-3-2, EN 61010
EEC guidelines: 89/336/EEC, 73/23/EEC

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Carrousel for
flame sterilizing
inoculation loops

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Flame time

Precise settings of the required flame sterilization time

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Cooling period

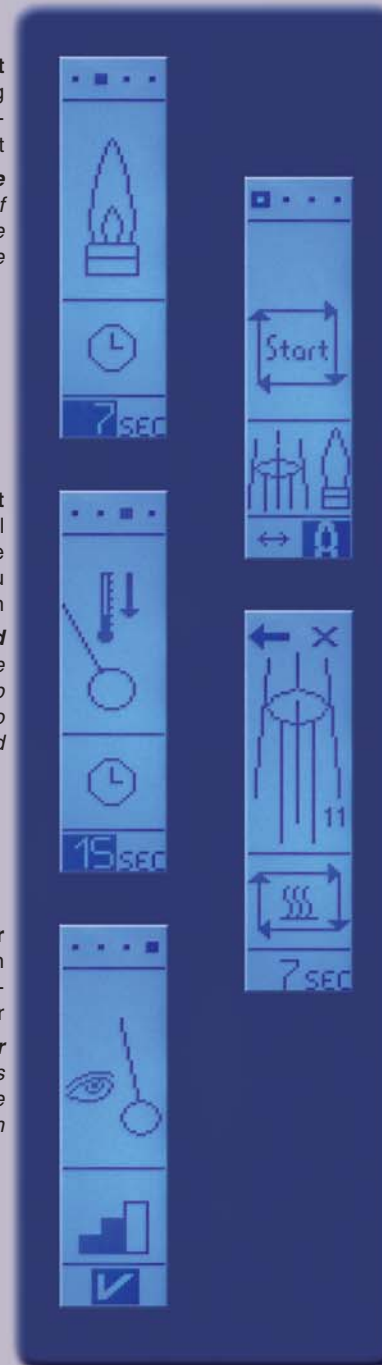
The cool-down time can be set for the inoculation loop material used - exact to the second

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Sterilization monitor

An intelligent sensor monitors the safe sterilization. Three setting options /off switch

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Start options

Start immediately or sterilize all inoculation loops in a full rotation

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Everything visible at a glance

Removal positions of cool inoculation loops or the remaining cool-down time, residual heat display and sterilization time

The fully automatic carousel for flame sterilizing inoculation loops

The inoculation loop carousel **Autoloop^{PRO}**, which is suitable for up to 4 inoculation loops, was developed to comply with current requirements in microbiological, cellbiological and biotechnological laboratories.

Modern Design

The stable housing enables comfortable and easy access to inoculation loops. Removal positions on both sides make the carousel equally suitable for right and left handers.

More Comfort

Keep all functions in view with the fully graphic display. The graphic menu control makes it easy to select all functions quickly. Flaming and cooling time can be adjusted to the second. During the flame sterilizing, the display shows the remaining flaming time. The **Autoloop^{PRO}** rotates and controls flaming automatically.

No unintentional use of hot inoculation loops: When flaming is completed, the display shows the remaining cooling time and the removal positions of cool inoculation loops.

All the suspended inoculation loops can be flamed as needed in one full rotation of the carousel.

More Safety

The nearly perpendicular position of the inoculation loop holder protects even the sleeve nut of the holder from possible contaminations. Additionally, an intelligent sensor of the **Autoloop^{PRO}** monitors safe sterilization.

Saves Energy, Time and Costs

Precise flame sterilization time extends the life of the inoculation loops and reduces gas consumption. Continuous working during the flaming and the cooling phases of the inoculation loops saves a great deal of time and makes it possible to work efficiently.

Robust

Fabricated entirely of stainless steel, anodized aluminium and a display, protected by heat resistant glass, the **Autoloop^{PRO}** can withstand extreme laboratory conditions.

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Easy to use

All parameters are easy to set; quick start button for immediate start

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Mount safely

Coordinated mounts and reliable optical detectors for inoculation loop holders

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Optimal sterilization

Sterilization monitor for safe sterilization of inoculation loops

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